HP PLUS TOWER NITROGEN GENERATOR





North America & South America contact:

Europe, Asia, Africa, & Australia/Oceania contact:


VICI DBS USA

tel: +1 713-263-6970 fax: +1 713-263-6971 web: www.vicidbs.com

VICI AG INTERNATIONAL

tel: Int + 41 41 925-6200 fax: Int + 41 41 925-6201 web: www.vicidbs.com

The VICI DBS° HP Plus Tower produces nitrogen by utilizing a combination of filtration and pressure swing adsorption (PSA) technology. Standard compressed air is filtered by high efficiency coalescing filters to remove all contaminants down to 5 micron. For ultra sensitive applications such as GC carrier and make-up gas, units also include the additional heated catalyst module to ensure hydrocarbons are removed to <0.1 ppm. The air then passes through two columns filled with a proprietary blended carbon molecular sieve (CMS) which adsorbs O2, CO2 and moisture. These are desorbed to the atmosphere during the pressure swing cycle leaving a supply of ultra pure nitrogen.

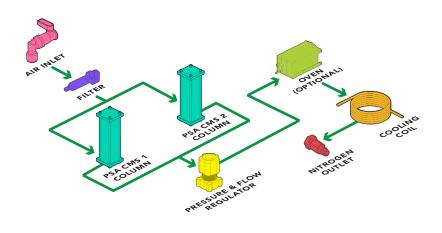



INCREASE EFFICIENCY

A constant gas supply with a guaranteed purity, eliminates interruptions of analyses to change cylinders and reduces the amount of instrument re-calibrations required.



RETURN ON INVESTMENT

Payback period can be as short as 6 to 12 months.



IMPROVE SAFETY

Nitrogen produced at low pressure and ambient temperature, removes the need for high pressure cylinders.



ENHANCE PERFORMANCE

Gas generators can be installed in the laboratory close to the instrument, eliminating the need for long gas lines from external cylinder supplies. A constant guaranteed high purity gas supply improves stability and ensures greater reproducibility of results.



FEATURES

Produces a continuous supply of high purity nitrogen | On-demand supply 24/7 | Flow rate: 200 to 4000 mL/min | Purity: up to +99.999% & <0.1 ppm THC | Pressure: up to 5 barg (75 psig) | Proprietary carbon molecular sieve technology | 2-year complete product warranty | Easy to install, operate and maintain



BENEFITS

Eliminates dangerous high pressure cylinders helping to keep your employees safer | Removes the logistics, inconvenience, downtime and costs of a cylinder system | Flow capacity to match your specific instrument demands | Ideal for all GC applications - stable baseline with increased sensitivity and repeatability | Meets and exceeds the requirements for the most demanding GC applications | Superior air purification with long life catalyst technology | Peace of mind | Improve your laboratory work flow and productivity



APPLICATIONS

GC APPLICATIONS

- GC carrier and make-up gas
- ECD
- ELSD

- TGA & DSC
- Incubators



| MODELS & SPECS | HP PLUS 500 | HP PLUS 750 | HP PLUS 1300 | HP PLUS 4000 | | | |
|---|---|-------------|---------------------|--------------|--|--|--|
| Flow mL/min | 500 | 750 | 1300 | 4000 | | | |
| Purity | +99. | 999% | +99.99% | +99% | | | |
| Hydrocarbon purity (measured as methane) | n/a | | | | | | |
| Dewpoint °C (°F) | -50 (-58) | | | | | | |
| Outlet pressure barg (psig) | up to 5 max (75) | | | | | | |
| Inlet pressure barg (psig) | 7 to 10 (100 to 160) | | | | | | |
| Actual inlet air requirement litres @ 8 barg (psig) | 11 (159.5) | 12 (174) | 16 (232.1) | 24 (348.1) | | | |
| Recommended compressor air inlet @ 8 barg (psig) | 22 (319.1) | 24 (348.1) | 32 (494.1) | 48 (696.2) | | | |
| Pressure drop barg (psig) | 1.5 (22) | | | | | | |
| Inlet air quality | Clean dry compressed air ISO8573-1:2010 Class 1.2.1 | | | | | | |
| Technology | Carbon molecular sieve | | | | | | |
| Warm up time minutes | 60 | | | | | | |
| LED indicators | Power on/off, system ready, errors | | | | | | |
| Electrical supply | 110-120V 60Hz / 220-240V 50 Hz | | | | | | |
| Power consumption watts | 12 | | | | | | |
| Noise level | Minimal | | | | | | |
| Dimensions mm (in) | 175W x 490H x 670D (6.9W x 19.3H x 26.4D) | | | | | | |
| Weight kg (lb) | 24.5 (54) | | | | | | |
| Shipping dimensions mm (in) | 770W x 410H x 590D (30.3W x 16.1H x 23.2D) | | | | | | |
| Shipping weight kg (lb) | 29.5 (65) | | | | | | |
| Operating temp °C (°F) | 15 to 35 (59 to 95) | | | | | | |
| Inlet connection | 1/4" Compression | | | | | | |
| Outlet connection | 1/8" Compression | | | | | | |
| Certification | CE, FCC | | | | | | |

| MODELS & SPECS | HP PLUS 200 HC | HP PLUS 500 HC | HP PLUS 750 HC | HP PLUS 1300 HC | HP PLUS 4000 HC | | |
|---|---|-------------------|-------------------|--------------------|--------------------|--|--|
| Flow mL/min | 200 | 500 | 750 | 1300 | 4000 | | |
| Purity | +99.999% +99.99% +99. | | | | | | |
| Hydrocarbon purity (measured as methane) | 0.1 ppm | | | | | | |
| Dewpoint °C (°F) | -50 (-58) | | | | | | |
| Outlet pressure barg (psig) | Up to 5 max (75) | | | | | | |
| Inlet pressure barg (psig) | 7 to 10 (100 to 160) | | | | | | |
| Actual inlet air requirement liters @ 8 barg (psig) | 11 (159.5) | 11 (159.5) | 12 (174) | 16 (232.1) | 24 (348.1) | | |
| Recommended compressor air inlet @ 8 barg (psig) | 22 (319.1) | 22 (319.1) | 24 (348.1) | 32 (494.1) | 48 (696.2) | | |
| Pressure drop barg (psig) | 1.5 (22) | | | | | | |
| Inlet air quality | Clean dry compressed air ISO8573-1:2010 Class 1.2.1 | | | | | | |
| Technology | Carbon molecular sieve | | | | | | |
| Warm up time minutes | 60 | | | | | | |
| LED indicators | Power on/off, system ready, errors | | | | | | |
| Electrical supply | 110-120V 60Hz / 220-240V 50 Hz | | | | | | |
| Power consumption watts | 270 | | | | | | |
| Noise level | Minimal | | | | | | |
| Dimensions mm (in) | 175W x 490H x 670D (6.9W x 19.3H x 26.4D) | | | | | | |
| Weight kg (lb) | 26.9 (59.3) | | | | | | |
| Shipping dimensions mm (in) | 770W x 410H x 590D (30.3W x 16.1H x 23.2D) | | | | | | |
| Shipping weight kg (lb) | 31.9 (70) | | | | | | |
| Operating temp °C (°F) | 15 to 35 (59 to 95) | | | | | | |
| Inlet connection | 1/4" Compression | | | | | | |
| Outlet connection | 1/8" Compression | | | | | | |
| Certification | CE, FCC | | | | | | |

ORDERING INFORMATION (for best service, please call to discuss your application before placing your order).

| HP PLUS 500 | | HP PLUS 4000 | | HP PLUS 750 HC | |
|---------------------|-----------|---------------------|-----------|------------------|-----------|
| DB-N2T-500-EU | 220V/50Hz | DB-N2T-4000-EU | 220V/50Hz | DB-N2T-750-O-EU | 220V/50Hz |
| DB-N2T-500-US | 115V/60Hz | DB-N2T-4000-US | 115V/60Hz | DB-N2T-750-O-US | 115V/60Hz |
| HP PLUS 750 | | HP PLUS 200 HC | | HP PLUS 1300 HC | |
| DB-N2T-750-EU | 220V/50Hz | DB-N2T-200-O-EU | 220V/50Hz | DB-N2T-1300-O-EU | 220V/50Hz |
| DB-N2T-750-US | 115V/60Hz | DB-N2T-200-O-US | 115V/60Hz | DB-N2T-1300-O-US | 115V/60Hz |
| HP PLUS 1300 | | HP PLUS 500 HC | | HP PLUS 4000 HC | |
| DB-N2T-1300-EU | 220V/50Hz | DB-N2T-500-O-EU | 220V/50Hz | DB-N2T-4000-O-EU | 220V/50Hz |
| DB-N2T-1300-US | 115V/60Hz | DB-N2T-500-O-US | 115V/60Hz | DB-N2T-4000-O-US | 115V/60Hz |